

Flexcell®'s Culture Plates Compatibility Guide

| When using the following culture plates | Required Loading Stations | Plates usable with the following systems if upgraded with the listed accessories |
|---|---|---|
| BioFlex® 6-well culture plates (page 19) | BioFlex® Loading Stations™ | <p>FX-5000T and FX-6000T require no further equipment</p> <p>FX-5000TT and FX-6000TT require BioFlex® Baseplate Kit or optional separately available BioFlex® Loading Stations™</p> <p>FX-5000C requires Tension FlexLink® with BioFlex® Baseplate and vacuum pump</p> |
| BioPress™ 6-well Compression culture plates (page 24) | Not required | <p>FX-5000T and FX-6000T require Compression FlexLink® with Compression Baseplate and compressor</p> <p>FX-5000TT and FX-6000TT require Compression FlexLink® with Compression Baseplate and compressor</p> <p>FX-5000C requires no further equipment</p> |
| HT BioFlex® 24-well culture plates (page 19) | 24-well Loading Stations™ | <p>FX-5000T and FX-6000T require 24-well HT Baseplate Kit</p> <p>FX-5000TT and FX-6000TT require 24-well HT Baseplate Kit</p> <p>FX-5000C requires Tension FlexLink®, vacuum pump and 24-well HT Baseplate Kit</p> |
| Tissue Train® 6-well culture plates (page 20) | Arctangle® Loading Stations™, Trough Loaders™ | <p>FX-5000T and FX-6000T require Tissue Train® Baseplate Kit or optional separately available Arctangle® Loading Stations™ and Trough Loaders™</p> <p>FX-5000TT and FX-6000TT require no further equipment</p> <p>FX-5000C requires Tension FlexLink®, vacuum pump and Tissue Train® Baseplate Kit, or Tension FlexLink®, vacuum pump and optional separately available Arctangle® Loading Stations™ and Trough Loaders™</p> |
| Tissue Train® Circular Foam 6-well culture plates (page 21) | BioFlex® Loading Stations™ | <p>FX-5000T and FX-6000T require no further equipment</p> <p>FX-5000TT and FX-6000TT require BioFlex® Baseplate Kit or optional separately available BioFlex® Loading Stations™</p> <p>FX-5000C requires Tension FlexLink® and vacuum pump</p> |
| Tissue Train® Trapezoidal 6-well culture plates (page 20) | Arctangle® Loading Stations™, Trapezoidal Trough Loaders™ | <p>FX-5000T and FX-6000T require Tissue Train® Baseplate Kit and Trapezoidal Trough Loaders™ or optional separately available Arctangle® Loading Stations™ and Trapezoidal Trough Loaders™</p> <p>FX-5000TT and FX-6000TT require Trapezoidal Trough Loaders™</p> <p>FX-5000C requires Tension FlexLink®, vacuum pump, Tissue Train® Baseplate Kit and Trapezoidal Trough Loaders™; or Tension FlexLink®, vacuum pump, optional separately available Arctangle® Loading Stations™ and Trapezoidal Trough Loaders™</p> |
| UniFlex® 6-well culture plates (page 21) | Arctangle® Loading Stations™ | <p>FX-5000T and FX-6000T require Tissue Train® Baseplate Kit or optional separately available Arctangle® Loading Stations™</p> <p>FX-5000TT and FX-6000TT require no further equipment</p> <p>FX-5000C requires Tension FlexLink®, vacuum pump and Tissue Train® Baseplate Kit; or Tension FlexLink®, vacuum pump and optional separately available Arctangle® Loading Stations™</p> |